

# EASYMIX

## TECHNICAL FEATURES AND INSTALLATION RECOMMENDATIONS

### A - TECHNICAL FEATURES

#### AESAP1068 EASYMIX BLENDER

#### AESAP10 EASYMIX BLENDER WITH GLASS DOOR



<b>Power input</b>	100 Watts maximum	
<b>Power supply</b>	115V +/- 10%, 50-60Hz or 230V +/- 10%, 50-60Hz, 70W, 0,5 / 0,3A	
<b>Calibre des fusibles</b>	2AT	
<b>Fuses rating</b>	From 5 to 40°C	
<b>Relative humidity</b>	From 50 to 80%	
<b>Atmospheric pressure</b>	From 700 HPa to 1100 HPa	
<b>Height</b>	Up to 2000 meters	
<b>Pollution level</b>	2	
<b>Installation type</b>	Type II according to IEC664 directives	
<b>Dimensions (WxHxD)</b>	Unit only (door closed): 260 x 265 x 400 mm	Packed 310 x 424 x 500 mm
<b>Weight</b>	Unit only: 20 Kg	
<b>Motor speed</b> (empty unit) (Value depends on power supply)	From 200 to 260 rpm	
<b>Number of paddle strokes / min</b> (empty unit) (Value depend on power supply)	From 400 to 520 strokes / min	
<b>Adjustable timer</b>	From 10 to 180 seconds or continuous	
<b>2x 16 character alphanumerical display ( with backlight)</b>		
<b>Optional wastage tray (AESMI0070)</b>		
<b>Dimensions (W x l x D mm) : 250 x 350 x 35</b>		

# EASYMIX

## TECHNICAL FEATURES AND INSTALLATION RECOMMENDATIONS

### B - INSTALLATION RECOMMENDATIONS

Easy MIX must be installed close to a wall socket.  
Vibrations are created while the unit is operating. It must be placed on a flat, stable and horizontal surface.  
It should be placed on a dry and clean workbench

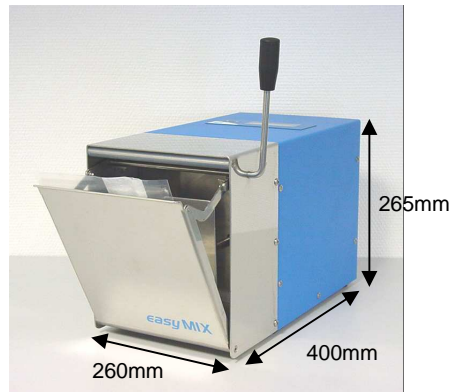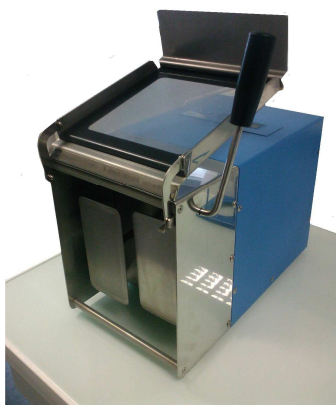

Only the **bent power cable** delivered with the unit must be used.  
AES Chemunex will not be responsible for any damage due to the use of a different cable.

Before plugging in the unit, make sure the voltage selector at the back is positioned according to the room mains power features.

When the unit is delivered, the cursor is placed to fit with a 230V +/-10% power supply.



### C - CONSOMMABLES ET ACCESORIES

CONSOMMABLES	
<b>AES400/50G</b>	Type 400 standard sterile irradiated bag (2500 units - 50 packs of 50 bags) <i>For optimal results use the bags recommended by AES CHEMUNEX</i>
<b>AES400P/50G</b>	Type 400 sterile bags with side filter (1500 units- 30 packets of 50 bags) <i>For optimal results use the bags recommended by AES CHEMUNEX</i>
<b>AES400FB/25G</b>	Full filter sterile bags (1500 units - 60 packs of 25 bags) <i>For optimal results use the bags recommended by AES CHEMUNEX</i>
ACCESSOIRES	
<b>AESAP0678</b>	Stand alone altuglass bag holder 1 bag
<b>AESAP0679</b>	Stand alone altuglass bag holder 10 bags
<b>DCPBF225B</b>	Bag closing clips (50 units)